



Annexin V-FITC Apoptosis Detection Kit

Manufactured by BioVision.

ALX-850-250-KI01: ~25 tests
ALX-850-250-KI02: ~100 tests
ALX-850-250-KI03: ~400 tests

(Version 01: December 1, 2004)

NORTH/SOUTH AMERICA

ENZO LIFE SCIENCES
INTERNATIONAL, INC.
5120 Butler Pike
Plymouth Meeting, PA 19462-1202
USA
Phone: 1-800-942-0430/(610) 941-0430
Fax: (610) 941-9252
E-mail: info-usa@enzolifesciences.com

GERMANY

ENZO LIFE SCIENCES GMBH
Marie-Curie-Strasse 8
DE-79539 Lörrach
Germany
Phone: +49/0 7621 5500 526
Fax: +49/0 7621 5500 527
E-mail: info-de@enzolifesciences.com

SWITZERLAND & REST OF EUROPE

ENZO LIFE SCIENCES AG
Industriestrasse 17, Postfach
CH-4415 Lausen
Switzerland
Phone: +41/0 61 926 89 89
Fax: +41/0 61 926 89 79
E-mail: info-ch@enzolifesciences.com

BENELUX

ENZO LIFE SCIENCES BVBA
Melkerijweg 3
BE-2240 Zandhoven
Belgium
Phone: +32/0 3 466 04 20
Fax: +32/0 3 466 04 29
E-mail: info-be@enzolifesciences.com

UK & IRELAND

ENZO LIFE SCIENCES (UK) LTD.
Palatine House
Matford Court
Exeter EX2 8NL
UK
Phone: 0845 601 1488 (UK customers)
+44/0 1392 825900 (from overseas)
Fax: +44/0 1392 825910
E-mail: info-uk@enzolifesciences.com

www.enzolifesciences.com

For laboratory use only. Not for human or diagnostic use.

I. Introduction:

The **Annexin V-FITC Apoptosis Detection Kit** is based on the observation that soon after initiating apoptosis, most cell types translocate the membrane phospholipid phosphatidylserine (PS) from the inner face of the plasma membrane to the cell surface. Once on the cell surface, PS can be easily detected by staining with a fluorescent conjugate of Annexin V, a protein that has a high affinity for PS. The one-step staining procedure takes only 10 minutes. In addition, the assay can be directly performed on live cells, without the need for fixation. Detection can be analyzed by flow cytometry or by fluorescence microscopy.

II. Kit Contents:

Components	850-250-KI01	850-250-KI02	850-250-KI03
	25 tests	100 tests	400 tests
Annexin V-FITC	125 µl	500 µl	2 ml
1X Binding Buffer	12.5 ml	50 ml	2 x 100 ml
Propidium Iodide (PI)	125 µl	500 µl	2 ml

III. Annexin V-FITC Assay Protocol:

A. Incubation of cells with Annexin V-FITC

1. Induce apoptosis by desired method.
2. Collect $1-5 \times 10^5$ cells by centrifugation.
3. Resuspend cells in 500 µl of 1x Binding Buffer.
4. Add 5 µl of Annexin V-FITC and 5µl of propidium iodide (PI, optional.)
5. Incubate at room temperature for 5 min in the dark.

Proceed to B or C below depending on method of analysis.

B. Quantification by Flow Cytometry

Analyze Annexin V-FITC binding by flow cytometry (Ex = 488 nm; Em = 530 nm) using FITC signal detector (usually FL1) and PI staining by the phycoerythrin emission signal detector (usually FL2).

For adherent cells, gently trypsinize and wash cells once with serum-containing media before incubation with Annexin V-FITC (A.3-5).

C. Detection by Fluorescence Microscopy

1. Place the cell suspension from Step A.5 on a glass slide. Cover the cells with a glass coverslip.

For analyzing adherent cells, grow cells directly on a coverslip. Following incubation (A.5), invert coverslip on a glass slide and visualize cells. The cells can also be washed and fixed in 2% formaldehyde before visualization. (Cells must be incubated with Annexin V-FITC before fixation since any cell membrane disruption can cause nonspecific binding of Annexin V to PS on the inner surface of the cell membrane).

2. Observe the cells under a fluorescence microscope using a dual filter set for FITC & rhodamine.

Cells that have bound Annexin V-FITC will show green staining in the plasma membrane. Cells which have lost membrane integrity will show red staining (PI) throughout the nucleus and a halo of green staining (FITC) on the cell surface (plasma membrane).

IV. Storage:

Store kit at +4°C. Avoid freeze/thaw cycles. Protect from light!